

ET 351C

Thermodynamics of the Refrigeration Circuit



- * **Compression refrigeration system for thermodynamic investigations**
- * **Indirectly heated evaporator and water-cooled condenser**
- * **Open compressor with pendulum bearing drive motor for torque measurement**

Technical Description

In this trainer, great care was taken to make the thermodynamic processes in the refrigeration system as transparent as possible. The capacities of compressor, evaporator and condenser can be measured. Pressure and temperature measuring points are provided at all the relevant locations to also allow for the pressure and heat losses in a refrigeration system to be investigated in detail.

The refrigeration circuit of ET 351C contains an open compressor with variable speed, a water-cooled condenser, a thermostatic expansion valve and an evaporator heated by a hot water circuit.

The compressor is driven via a pendulum bearing motor with frequency converter for speed adjustment. A force transducer permits the measuring of the drive torque. Using the speed the mechanical drive power of the compressor can thus be determined. The electrical heating power of the hot water circuit can be freely adjusted and is displayed. The condenser capacity is measured via the cooling water flow.

The measured values can be read on digital displays. At the same time, the measured values can also be transmitted directly to a PC via USB. The data acquisition software is included. The software allows for the process to be represented in the log p-h diagram and displays the key characteristic variables of the process, e.g. compressor pressure ratio and coefficient of performance.

The well-structured instructional material sets out the fundamentals and provides a step-by-step guide through the experiments.

Learning Objectives / Experiments


- cyclic process in the log p-h diagram
- comparison of the real cyclic process and the ideal cyclic process
- energy balances at the evaporator and condenser
- calculation of the motor power via speed and torque
- determination of losses
- calculation of the coefficient of performance
- operating behaviour under load
- non-steady-state operating behaviour

ET 351C


Thermodynamics of the Refrigeration Circuit



1 expansion valve, 2 evaporator, 3 refrigerant flow meter, 4 pressure switch, 5 process schematic, 6 receiver, 7 hot water circuit of the evaporator, 8 drive motor, 9 compressor, 10 cooling water flow meter, 11 condenser, 12 displays and controls

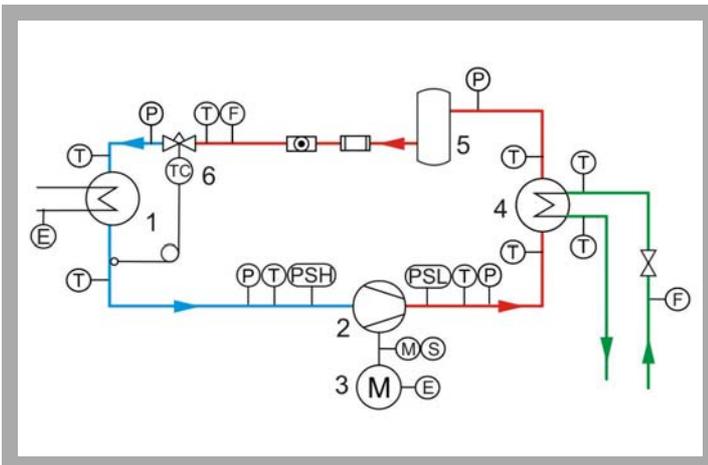

1 evaporator, 2 compressor, 3 drive motor, 4 condenser, 5 receiver, 6 expansion valve; P pressure, T temperature, M torque, S speed, E electrical power; PSL, PSH pressure switch; blue: low pressure, red: high pressure, green: cooling water

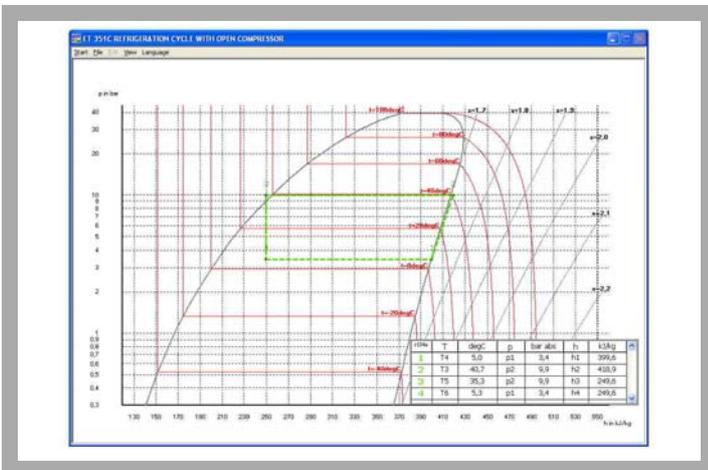

Software screenshot: log p-h diagram

Specification

- [1] thermodynamic investigation of a refrigeration circuit
- [2] refrigeration circuit with open compressor, water-cooled condenser, thermostatic expansion valve and indirectly heated evaporator
- [3] compressor drive with speed-controlled motor via V-belt
- [4] motor on pendulum bearing for torque measurement
- [5] tube evaporator with hot water circuit as cooling load
- [6] water-cooled coaxial coil heat exchanger as condenser
- [7] displays for temperature, pressure, flow rate, speed, torque and power at the equipment
- [8] GUNT software for data acquisition via USB under Windows Vista or Windows 7
- [9] refrigerant R134a, CFC-free

Technical Data

- Open compressor
 - refrigeration capacity: approx. 550W (at a speed of 500min⁻¹ and -10/20°C)
 - motor: 550W, 1.400min⁻¹
- Heater: 1x 1.000W
- Condenser
 - capacity: 1.300W
- Measuring ranges
 - temperature: 9x -30...100°C, 1x 0...100°C
 - pressure: 1x -1...9bar, 1x -1...24bar, 2x -1...15bar
 - torque (compressor): 0...10Nm
 - speed (compressor): 0...2.500min⁻¹
 - power consumption (compressor): 0...1.125W
 - power (heater): 0...1.125W
 - flow rate (water): 5...70g/s
 - flow rate (refrigerant): 0...0,5L/min

Dimensions and Weight

LxWxH: 1.520x790x1.760mm
Weight: approx. 120kg

Required for Operation

230V, 50/60Hz, 1 phase
Water connection

Scope of Delivery

- 1 trainer
- 1 set of hoses, 1 manual pump
- 1 GUNT software CD + USB cable
- 1 set of instructional material

Order Details

061.351C0 ET 351C Thermodynamics of the Refrigeration Circuit